## Aim of the experiment

Systematic extraction, isolation, purification and identification of **non-metallic poison** by various analytical techniques.

## Requirements

- a) **Test sample:** Food curry (divided in four parts) spiked with sodium salts containing chloride, thiosulfate, nitrate and nitrite, respectively.
- b) **Materials:** Beakers, test-tubes, test-tube holder, test-tube, stands, Dropper, Pipette, Mortar-pestle, filter paper sheet.
- c) Instruments: Hot plate
- d) **Chemicals:** Hydrochloric acid, ferric chloride, sulfanilic acid, alpha-naphthylamine, zinc dust.

# **Theory and Principle**

- a. Non-metals are the elements which form negative ions by accepting or gaining electrons.
- b. The qualitative analysis of anions involves the detection and identification of radicals present in inorganic salts
- c. Chlorine is a halogen, an element belonging to Group 7 of periodic table (Chloride: halide ion). Halide ions are detected using silver nitrate solutions.
- d. Sodium thiosulfate with ferric chloride reacts to form ferrous thiosulfate (dark violet).
- e. 1-napthylamine and sulfanilic acid is used for detection of nitrate and nitrite where reductase reduces nitrate to nitrite and reacts with sulfanilic acid to form diazonium salt. This salt on reaction with 1-napthylamine forms an azo dye.

#### **Procedure**

# **Extraction of poison:**

- a) Add 50 ml of hydrochloric acid (for digestion of sample) to the sample matrix in beaker along with equal amount of water.
- b) Keep the beaker(s) on hot plate to heat for 20-30 minutes (or till digestion was completed) at 110 degrees celsius.
- c) During heating HCL can be added if loss of volume is observed.
- d) The beaker(s) is then left to cool. Further, digested sample is filtered and filtrate can be used for anion analysis.

# **Reagent Preparation:**

- a) Preparation of alpha-naphthylamine solution: 0.5 g α-naphthylamine to be dissolved in 5N acetic acid by gently heating and made up to 100 ml. 5N acetic acid solution can be prepared by adding distilled water to 28.7 ml glacial acetic acid (17.4 N) to give a final volume of 100 ml. OR
  - Prepared by boiling 0.3 gm of 1-napthylamine with 70 ml of water, filtering or decanting from the small residue and mixing with 30 ml of glacial acetic acid.
- b) Preparation of sulfanilic acid solution: Prepared by dissolving 1 gm of sulfanilic acid in 30% warm acetic acid and volume made up to 100 ml.

**Sample Preparation:** Acidified solution of extract was prepared by adding few drops of acetic acid to be used for nitrate and nitrite test.

## **Detection of anions:**

Test	Procedure	Observation	Inference		
sample					
S1	A few drops acidified solution of the	Red colour will obtain.	Indicates presence of nitrate.		
	extracted sample was taken in a test-tube				
	+ 1 drop of sulfanilic acid				
	+ 1 drop of α-napthylamine				
	+ few mg of zinc dust was added				
S2	A few drops acidified solution of the	Red colour will obtain.	Indicates presence		
	extracted sample was taken in a test-tube		of nitrite		
	+ 1 drop of sulfanilic acid				
	+ 1 drop of α-napthylamine				
S3	2 ml extracted sample was taken in a	Dark Violet Colour	Indicates presence		
	test-tube	appears.	of thiosulphate.		
	+ Few drops of ferric chloride solution is				
	added.				
	On standing colour disappears.				
S4	Extracted sample was taken in a test-tube	White precipitate forms.	Indicates presence		
	+ Silver nitrate solution (1:1)		of chloride		
	Precipitate is soluble in dilute nitric acid but				
	soluble in dilute ammonia / potassium				
	cyanide / sodium thiosulphate				

# **Conclusion:**

On	the	basis	of	observation,	it	can	be	concluded	that	the	test	sample
respectively.												