Date:

Aim of the Experiment:

SAMPLING AND ISOLATION OF PROBIOTIC MICROORGANISMS

Introduction

Probiotics used by humans are usually isolated from different environments that include dairy and non-dairy sources. However, probiotics that are isolated from human or animal intestines have certain characteristics that differ from those isolated from dairy products. Isolation of *Lactobacillus* can be done from the milk sample by using selective media. *Lactobacillus* sp. was isolated from the milk samples by streak plate technique on MRS agar plates.

Materials

- 1. Samples for isolation (human breast milk, faecal samples, dairy products and fermented foods)
- 2. Inoculation needle
- 3. Incubator (set at 37°C)
- 4. Sterile test tubes (10 ml capacity)
- 5. Test tube stand
- 6. Sterile pipettes
- 7. Petri dishes

Procedure

A. Sampling

Collection and processing of faecal samples obtained from infants and adult

- 1. Faecal samples can be obtained from healthier adult volunteers and infants.
- 2. Samples were collected in a 5 mL sterile screw cap vial containing 0.1% peptone water.
- 3. After, bringing samples to laboratory nearly 1g of samples was transferred to 10 mL LAMBAV broth for enrichment containing 0.5 mg/mL LA.

Collection and processing of dahi and lassi samples

- 1. Samples of dahi and lassi were obtained from commercial dairy shops of nearby localities and also from traditional milk vendors.
- 2. Nearly, 1g of dahi and 1 mL of lassi was added to 0.1%(w/v) peptone water and mix by doing vigorous vortexing.
- 3. One millilitre of homogenized samples was added to 9 mL of LAMVAB broth for enrichment containing 0.5 mg/mL LA.

Collection and processing of mother breast milk samples

- 1. Mother's breast milk samples were collected with the help of trained nurse staff of the hospital in sterile 5 mL vials and transported immediately to laboratory in an ice box container.
- 2. Then 1 mL sample was enriched in 9.0 mL of LAMBAV broth containing 0.5 mg/mL LA.

Processing of Standard cultures

1. The received ampoules were aseptically broken, transferred to litmus chalk milk and then incubated at 37°C for 24-48 h. Next, these were transferred to MRS broth, incubated and checked for purity before employed in experiments.

B. Enumeration Procedure

- 1. Each sample was enriched at 37°C in modified MRS containing 0.5 mg/mL LA under micro-aerophilic conditions for 24 h.
- 2. After enrichment, 1.0 mL of sample was added to 9.0 mL of 0.1% (w/v) peptone water, vortexed and then serially diluted up to 10^{-4} - 10^{-10} dilutions.
- 3. To obtain colonies, 0.1 mL aliquot from appropriate dilutions was pour plated on MRS agar and further incubated at 37°C for 48 h.
- 4. Typical pin pointed yellow-colored colonies were randomly selected and transferred to MRS broth.
- 5. Further, cultures were checked for Gram and negative staining and thereafter, examined microscopically for preliminary purification and morphology.

Observations

A) Colony Morphology

Interpretation

B) Total microbial Load

Interpretation